LAB-V

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LAB-V

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Thank you for choosing the professional high precision scale LAB-V. All the features of this product were designed to the state of the art and are optimised for simple and straightforward use. If you any queries or problems not indicated in the operating manual, please contact our service.

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*The scale is equipped with a level indicator on the left bottom side of the front panel and four adjustable leveling feet.
Adjust the leveling feet until the bubble appears in the center circle of the indicator*

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ENGLISH

INTENDED USE

The LAB-V scale is used for weighing and counting products. This scale is intended for internal operations. The scale complies with the type described in the EC type Approval Certificate as well as the applicable requirement of EC Directives 89/336/EEC and 73/23/EEC.

SAFETY INSTRUCTIONS

Before using the scale, please read with care the information given in the operating manual. It contains important instructions for installation, proper use and maintenance of the appliance.

B3C SAS shall not liable for damages arising out of failure to heed the following instructions: When using electrical components under increased safety requirement, always comply with the appropriate regulations. Never perform work on the appliance while the power is switched on.

Improper installation will render the warranty null and void. Ensure the voltage marked on the power supply unit matches your mains power supply. The device meets the requirement for electromagnetic compatibility. Do not exceed the maximum values specified in the applicable standards.

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ENGLISH

CLEANING

It's sufficient to clean the scale with a moist cloth and a regular cleaning agent. Do not use any scourers.

MAINTENANCE

The scale does not require any routine maintenance. However, we recommend to check the scale's precision at regular intervals.

**NOTE**

Interference suppression of the unit corresponds to the valid EU directive 89/336/EEC. The display value may be influenced by extreme electrostatic and electro-magnetic interference, e.g. operation of a radio unit or a mobile phone in the immediate vicinity of the unit. When the electrostatic interference is no longer present, the unit can again be used for its intended purpose. You may need to switch it on again. If there is permanent electrostatic interference, we recommend to connect the platform with the ground. The scale is a measuring instrument. Air currents, vibrations, rapid changes in temperature and direct sunlight may have an effect on the accuracy of the scale.

WARRANTY - LIABILITY

This equipment is a measuring device.

No responsibility shall be accepted for damage caused through any of the following reasons: unsuitable or improper storage or use, incorrect or commissioning by the owner or third parties, natural wear, changes or modifications, incorrect or negligent handling, overuse, chemical, electrochemical or electrical interference. If operating, climatic or any other influences lead to a major change in conditions or material quality, the warranty for perfect until functioning shall be rendered null and void. Always keep the original packaging in case you have to return the scale.

NEDERLANDS**REINIGING**

It's sufficient to clean the scale with a moist cloth and a regular cleaning agent. Do not use any scourers.

ONDERHOUD

The scale does not require any routine maintenance. However, we recommend to check the scale's precision at regular intervals.

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▼ 2 - ECRAN / DISPLAY / DISPLAYTEKENS

15kg - 30kg - 50kg

OK Lecture stable / Reading is stable / Stabiele aflezing

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Unable Opération impossible / Unable requested operation / Operatie onmogelijk

HHHHH Surcharge / Weight exceed scale capacity / Overbelasten

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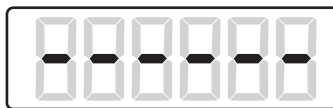
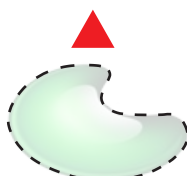
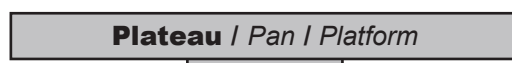
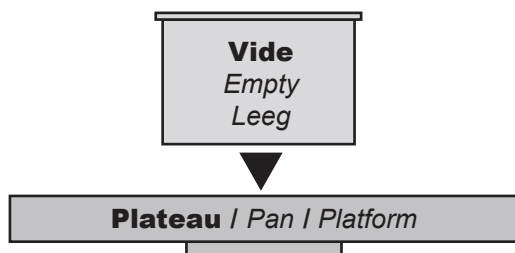
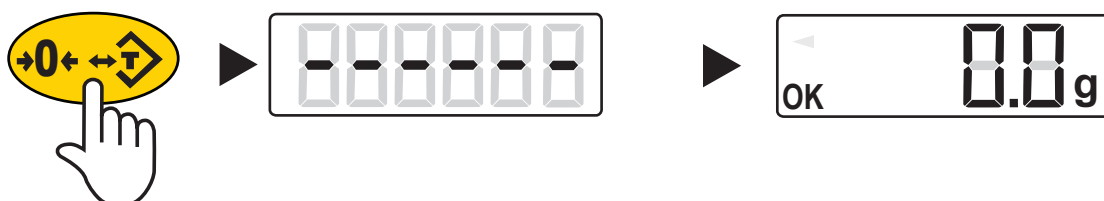
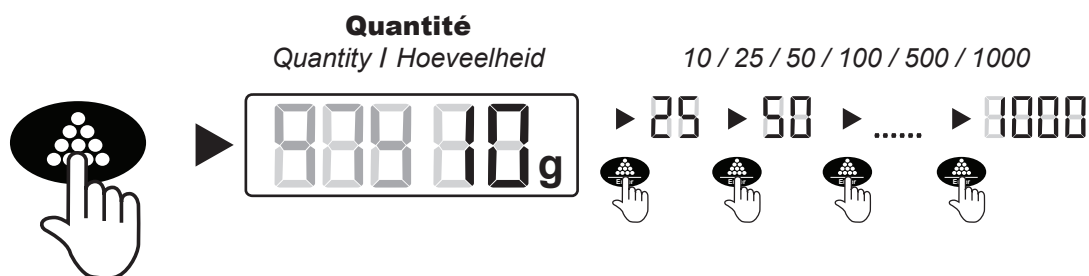
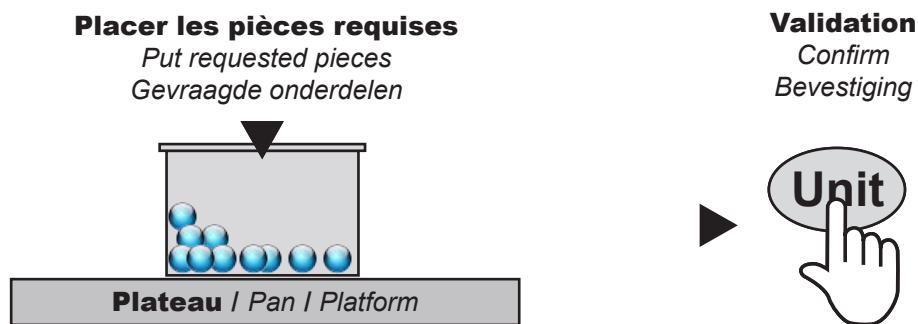
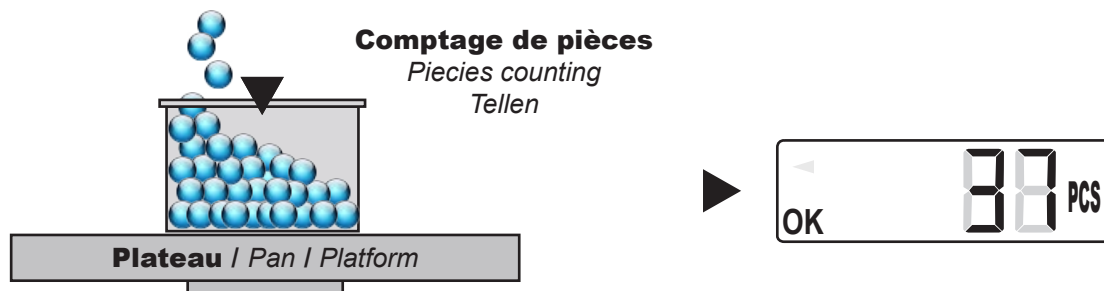
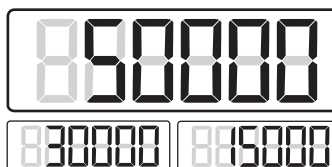
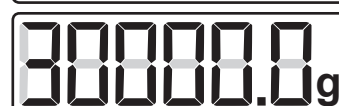
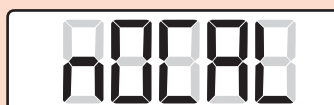
To calibrate at one half of the full capacity

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
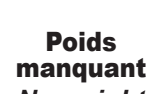
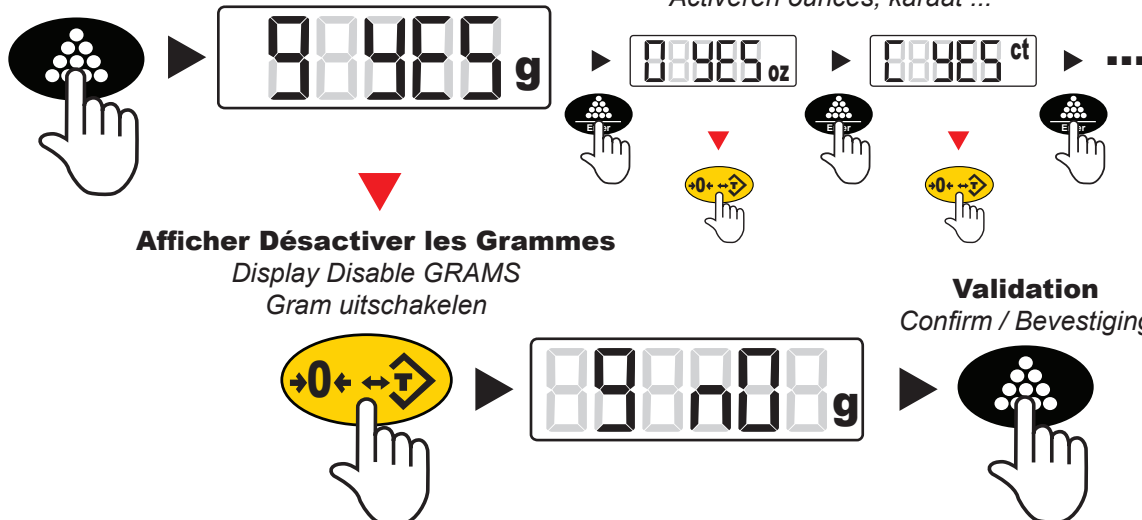
Return to weighing after the last unit

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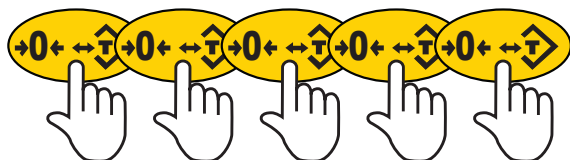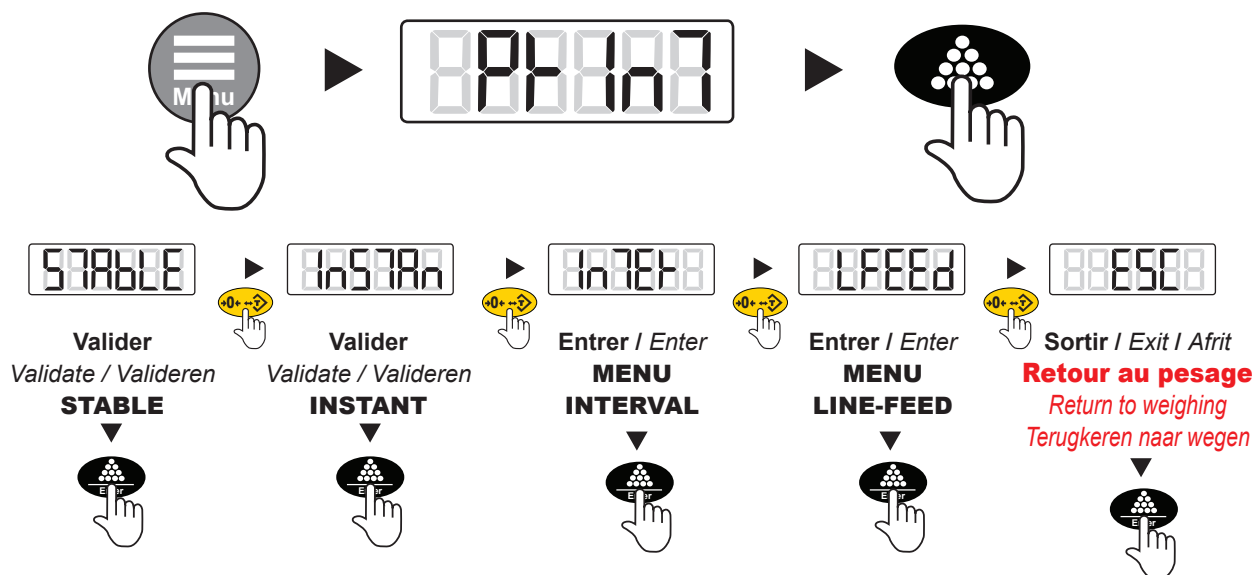
Once after weight change on standstill

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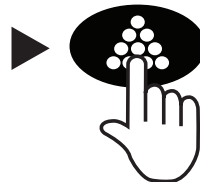
Continuous transmission (transmission rate adjustable)

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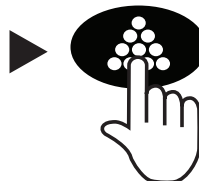
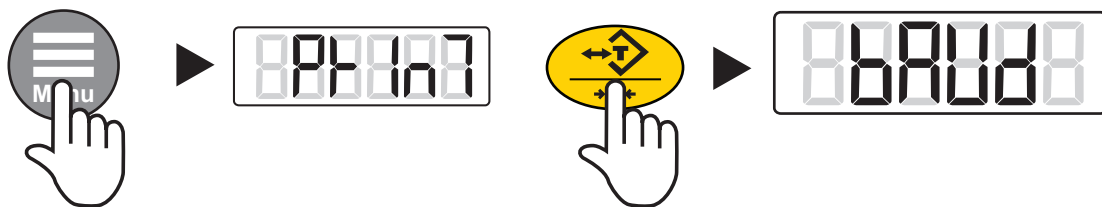
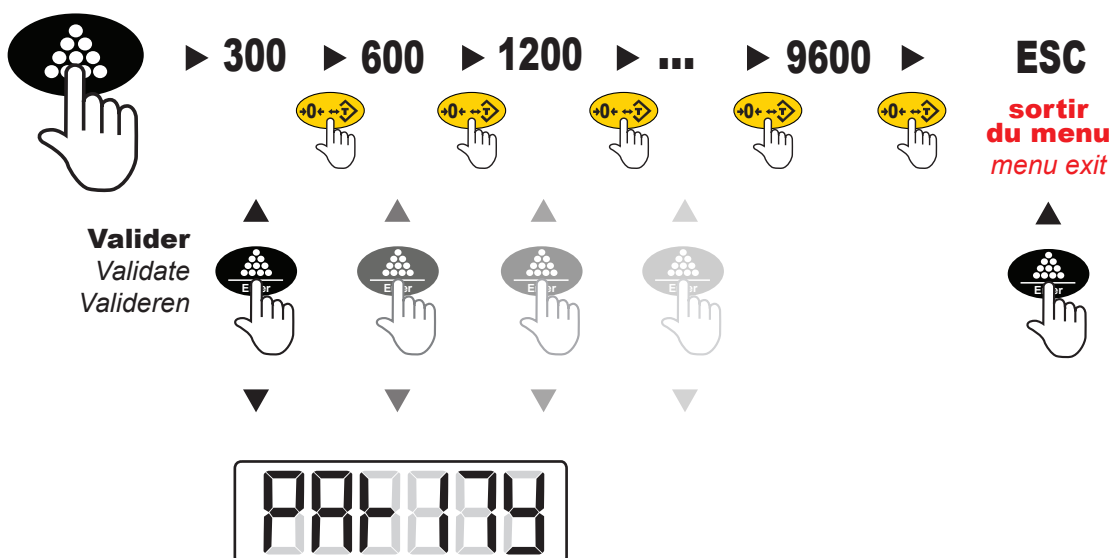
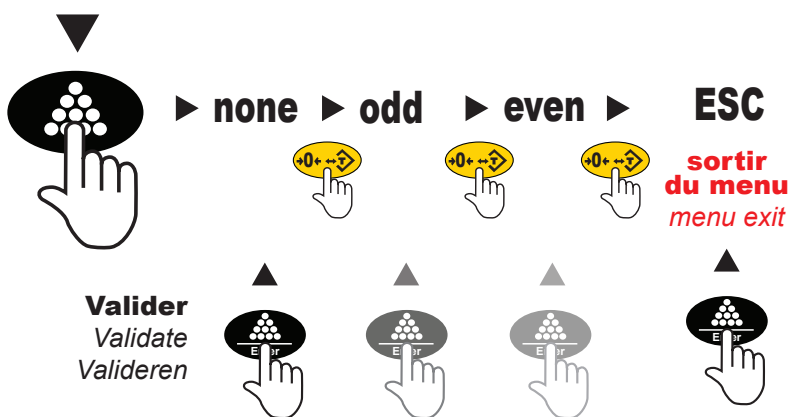
ENGLISH

Your balance can communicate with almost any RS232 device, the build-in interface does not include the complete protocol. Only the transmit and receive lines of the standard interface are used. This should not present any interfacing problems in most applications.

The data format is

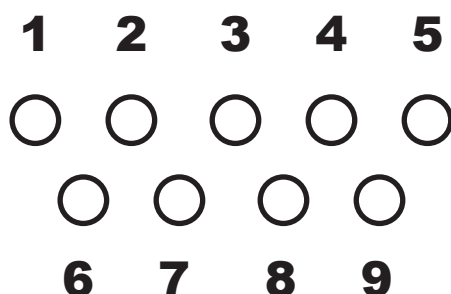
- 1 - start bit
- 8 - data bit include parity
- 1 - stop bit
- 10 - bit per frame (framing errors ignored)

Note: the balance will transmit using the parity selected.

However it does not check the parity it receives.

Use an RS232 cable to connect the external device to the balance, or construct one following the instructions below.

Connect a high quality, shielded cable with a DB9S (D-subminiature 9 pin female connector) using the following printout:

pin	description
2	TXD scale transmit data
3	RXD scale receives data
5	RRD signal ground

Note: “handshake” signals, such as “clear to send” (CES) are not used. The peripheral must have a minimum buffer (15 characters).

▼ NOTES...

b3c
pesage

Sérénité

www.b3c-pesage.fr

PESAGE - WEIGHING - WEGING

les justes valeurs, the true values, de echte waarden...

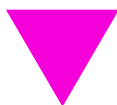
BOULANGERIE PATISSERIE

*Bakery - Pastry
Bakkerij - Gebak*

CUISINE PROFESSIONNELLE

*Professional kitchen
Professionele
keuken*

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IT - Bilance elettronica LAB-V

ND - Electronische weegschaal Reeks LAB-V

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The indicated balances respect the following European Standards

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